


# Clemex ST Mini 2.0 CCD Color Camera

USB 2.0 high speed camera installation sheet

**Note:** *The Clemex software must be installed before installing the camera.*




1. Please ensure that you have a **High Speed USB 2.0** port on your PC.
2. Right-click on your Desktop to access the **Screen Resolution** Window and check that your screen resolution is set at a minimum of 1920 x 1080.

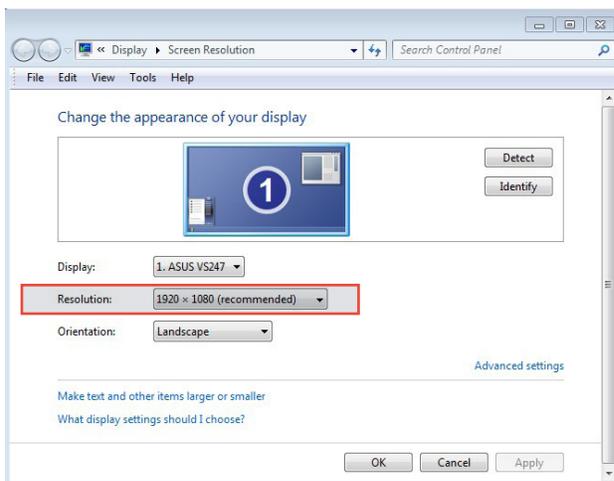



3. Plug the Clemex ST Mini 2.0 CCD Color camera in any available USB 2.0 port.

**CAUTION**  
*If you connect your camera to a non USB 2.0 USB port  
your camera may not work.*

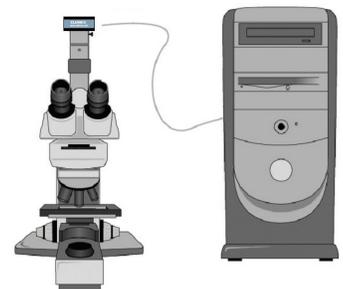

# Clemex ST Mini 2.0 CCD Color Camera


- The computer installs the appropriate device driver software and tells you when it is finished.

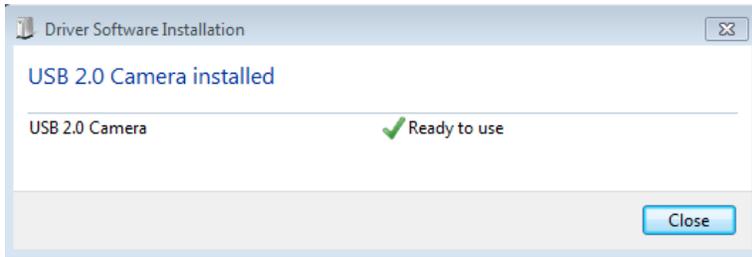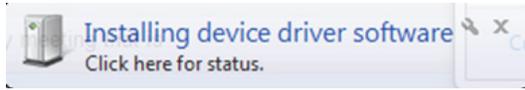

- To synchronise the camera with the Clemex software, open the **Video Input Setup** program which is found in the **Clemex Image Analysis** folder on your desktop.



- Click **Auto-Install Camera**. The camera details appear. Click to close the window.

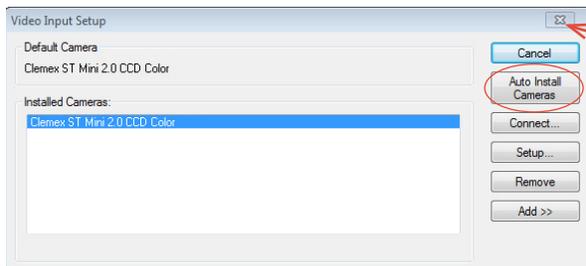

- Run your **Clemex Software** by double-clicking on the icon in the Clemex Software folder on your desktop.



Clemex Software